

Mouse Mitochondrial DNA Copy Number Assay Kit

(44 reactions)

Catalog Number: MCN 3

Store at -20°C.

FOR RESEARCH USE ONLY



Introduction: This DNA analysis kit is for the determination of mouse mitochondrial DNA copy number, *in vivo* and *in vitro*, by the comparison of mitochondrial (mt) and nuclear (n) DNA measured by real-time PCR.

Kit Contents:

- 96 well PCR plate
- rtPCR reaction mix.
- Validated primers (10 µM) to quantify mitochondrial DNA (mtDNA).
- Validated primers (10 µM) to quantify nuclear DNA (nDNA).
- Positive control [2.5 ng/ µl] (isolated total DNA from liver of B6 mouse).

Not Included in Kit:

- DNA isolation Kit
- Nuclease-free water
- PCR Tubes and Caps

Thermal cycler program:

- Preprogram PCR machine for this profile:
 - a. 95°C, 10 min
(40 Cycles)
 - b. 95°C, 15 sec
 - c. 60°C, 60 sec

Real time PCR procedure: The following procedure is for each 20 µL reaction. Increase all amounts proportionally according to the total number of tubes.

- Per PCR tube (20 µL Rx), mix the following:
 - a. 1 µL forward primer
 - b. 1 µL reverse primer
 - c. 8 µL sample contain genomic DNA/ 8 µL of positive control
 - d. 10 µL rtPCR reaction mix

Recommended concentration: Between 0.3 to 5.0 ng/µL


How to Calculate Mitochondrial Copy Number:

$$\Delta Ct1 = Ct (\text{mitochondria-control}) - Ct (\text{nucleus-control})$$

$$\Delta Ct2 = Ct (\text{mitochondrial-experimental}) - Ct (\text{nucleus-experimental})$$


$\Delta\Delta Ct = \text{Sample } \Delta Ct - \text{average } \Delta Ct \text{ control}$
 mtDNA fold change = $2^{-\Delta\Delta Ct}$

Total DNA isolated from mouse liver

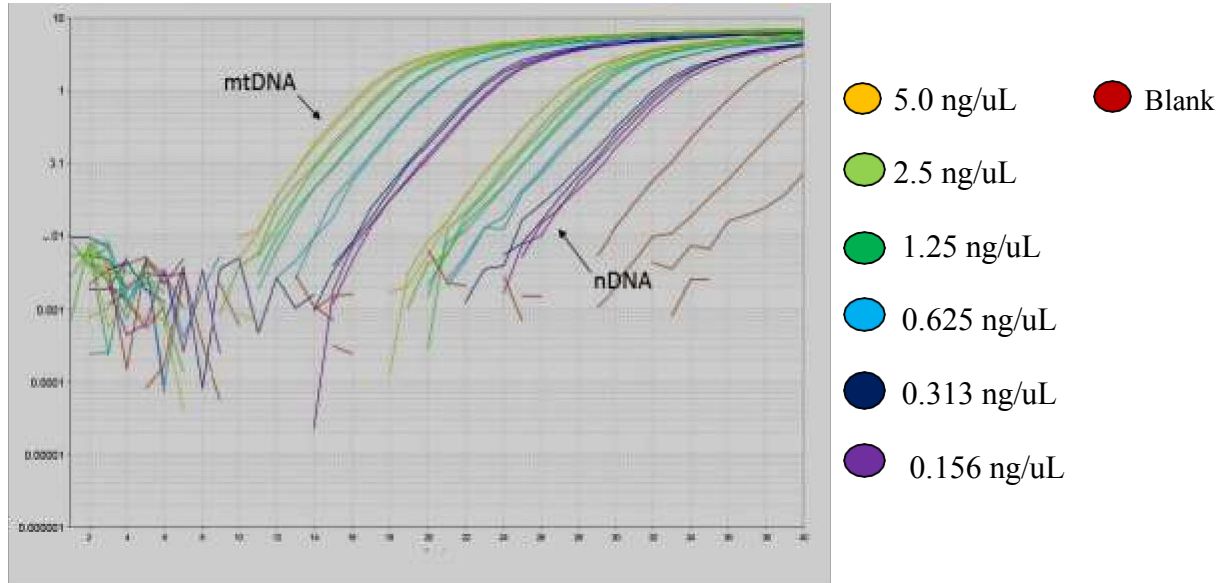

Suggested assay plate layout: n = nucleus; mt = mitochondria; BLK = blank

	1	2	3	4	5	6	7	8	9	10	11	12
A	nBLK	nS3	nS7	nS11	nS15	nS19	mtBLK	mtS3	mtS7	mtS11	mtS15	mtS19
B	nBLK	nS3	nS7	nS11	nS15	nS19	mtBLK	mtS3	mtS7	mtS11	mtS15	mtS19
C	nPC	nS4	nS8	nS12	nS16	nS20	mtPC	mtS4	mtS8	mtS12	mtS16	mtS20
D	nPC	nS4	nS8	nS12	nS16	nS20	mtPC	mtS4	mtS8	mtS12	mtS16	mtS20
E	nS1	nS5	nS9	nS13	nS17	nS21	mtS1	mtS5	mtS9	mtS13	mtS17	mtS21
F	nS1	nS5	nS9	nS13	nS17	nS21	mtS1	mtS5	mtS9	mtS13	mtS17	mtS21
G	nS2	nS6	nS10	nS14	nS18	nS22	mtS2	mtS6	mtS10	mtS14	mtS18	mtS22
H	nS2	nS6	nS10	nS14	nS18	nS22	mtS2	mtS6	mtS10	mtS14	mtS18	mtS22

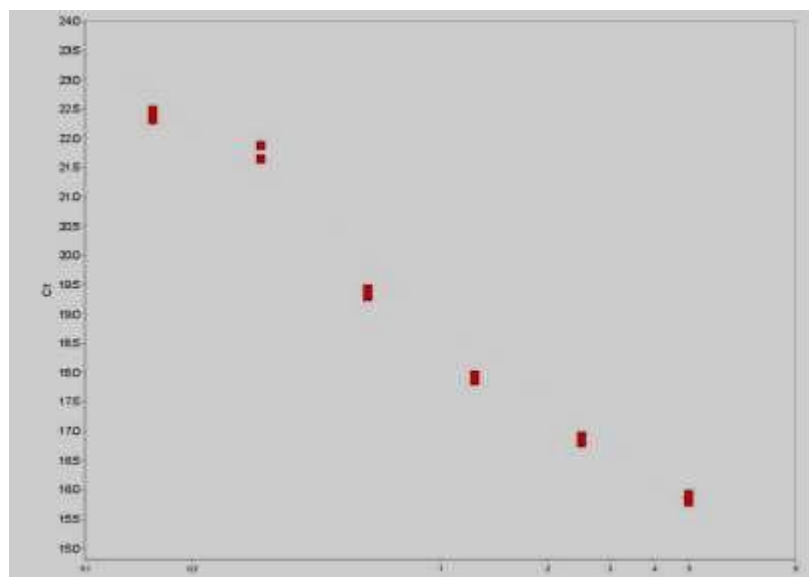Plot of C_T versus DNA concentration

References

1. Yi CX, Walter M, Gao Y, Pitra S, Legutko B, Kälın S, Layritz C, García-Cáceres C, Bielohuby M, Bidlingmaier M, Woods SC, Ghanem A, Conzelmann KK, Stern JE, Jastroch M, Tschöp MH. TNF α drives mitochondrial stress in POMC neurons in obesity. *Nat Commun*. 2017 May 10;8:15143. doi: 10.1038/ncomms15143. PMID: 28489068.
<https://www.ncbi.nlm.nih.gov/pmc/articles/PMC5436136/pdf/ncomms15143.pdf>
2. Marlein CR, Zaitseva L, Piddock RE, Robinson SD, Edwards DR, Shafat MS, Zhou Z, Lawes M, Bowles KM, Rushworth SA. NADPH oxidase-2 derived superoxide drives mitochondrial transfer from bone marrow stromal cells to leukemic blasts. *Blood*. 2017 Oct 5;130(14):1649-1660. doi: 10.1182/blood-2017-03-772939. PMID: 28733324.
<https://ashpublications.org/blood/article/130/14/1649/107660/NADPH-oxidase-2-derived-superoxide-drives>
3. Jessop F, Schwarz B, Heitmann E, Buntyn R, Wehrly T, Bosio CM. Temporal Manipulation of Mitochondrial Function by Virulent *Francisella tularensis* To Limit Inflammation and Control Cell Death. *Infect Immun*. 2018 Jul 23;86(8):e00044-18. doi: 10.1128/IAI.00044-18. PMID: 29760217.
<https://www.ncbi.nlm.nih.gov/pmc/articles/PMC6056872/pdf/e00044-18.pdf>
4. Singh SP, McClung JA, Bellner L, Cao J, Waldman M, Schragenheim J, Arad M, Hochhauser E, Falck JR, Weingarten JA, Peterson SJ, Abraham NG. CYP-450 Epoxygenase Derived Epoxyeicosatrienoic Acid Contribute To Reversal of Heart Failure in Obesity-Induced Diabetic Cardiomyopathy *via* PGC-1 α Activation. *Cardiovasc Pharm Open Access*. 2018;7(1):233. doi: 10.4172/2329-6607.1000233. PMID: 29707604.
<https://www.ncbi.nlm.nih.gov/pmc/articles/PMC5922773/pdf/nihms954607.pdf>

5. Basheer WA, Fu Y, Shimura D, Xiao S, Agvanian S, Hernandez DM, Hitzeman TC, Hong T, Shaw RM. Stress response protein GJA1-20k promotes mitochondrial biogenesis, metabolic quiescence, and cardioprotection against ischemia/reperfusion injury. *JCI Insight*. 2018 Oct 18;3(20):e121900. doi: 10.1172/jci.insight.121900. PMID: 30333316.
<https://www.ncbi.nlm.nih.gov/pmc/articles/PMC6237442/pdf/jciinsight-3-121900.pdf>
6. Helmer RA, Kaur G, Smith LA, Chilton BS. Helicase-like transcription factor (Hltf) gene-deletion promotes oxidative phosphorylation (OXPHOS) in colorectal tumors of AOM/DSS-treated mice. *PLoS One*. 2019 Aug 28;14(8):e0221751. doi: 10.1371/journal.pone.0221751. PMID: 31461471.
<https://www.ncbi.nlm.nih.gov/pmc/articles/PMC6713344/pdf/pone.0221751.pdf>
7. Akhtar S, Siragy HM. Pro-renin receptor suppresses mitochondrial biogenesis and function via AMPK/SIRT-1/ PGC-1 α pathway in diabetic kidney. *PLoS One*. 2019 Dec 4;14(12):e0225728. doi: 10.1371/journal.pone.0225728. PMID: 31800607.
<https://www.ncbi.nlm.nih.gov/pmc/articles/PMC6892478/pdf/pone.0225728.pdf>
8. Hou X, Li Z, Higashi Y, Delafontaine P, Sukhanov S. Insulin-Like Growth Factor I Prevents Cellular Aging via Activation of Mitophagy. *J Aging Res*. 2020 Aug 1;2020:4939310. doi: 10.1155/2020/4939310. PMID: 32802505.
<https://www.ncbi.nlm.nih.gov/pmc/articles/PMC7416301/pdf/JAR2020-4939310.pdf>
9. Tyrrell DJ, Blin MG, Song J, Wood SC, Goldstein DR. Aging Impairs Mitochondrial Function and Mitophagy and Elevates Interleukin 6 Within the Cerebral Vasculature. *J Am Heart Assoc*. 2020 Dec;9(23):e017820. doi: 10.1161/JAHA.120.017820. Epub 2020 Nov 23. PMID: 33225820.
<https://www.ncbi.nlm.nih.gov/pmc/articles/PMC7763766/pdf/JAH3-9-e017820.pdf>
10. Shimura D, Nuebel E, Baum R, Valdez SE, Xiao S, Warren JS, Palatinus JA, Hong T, Rutter J, Shaw RM. Protective mitochondrial fission induced by stress-responsive protein GJA1-20k. *Elife*. 2021 Oct 5;10:e69207. doi: 10.7554/eLife.69207. PMID: 34608863.
<https://www.ncbi.nlm.nih.gov/pmc/articles/PMC8492060/pdf/elife-69207.pdf>
11. Yuan Q, Lv Y, Ding H, Ke Q, Shi C, Luo J, Jiang L, Yang J, Zhou Y. CPT1 α maintains phenotype of tubules via mitochondrial respiration during kidney injury and repair. *Cell Death Dis*. 2021 Aug 14;12(8):792. doi: 10.1038/s41419-021-04085-w. PMID: 34392303.
https://www.ncbi.nlm.nih.gov/pmc/articles/PMC8364553/pdf/41419_2021_Article_4085.pdf